

BGO

Bismuth Germanate

Scintillation Material

Bismuth Germanate (BGO) is a high Z, high density scintillation material with chemical composition $\text{Bi}_4\text{Ge}_3\text{O}_{12}$. Due to the high atomic number of bismuth (83) and its high density, BGO is a very efficient γ -ray absorber.

BGO is a relatively hard, rugged, non-hygroscopic crystal which does not cleave. The material does not show any significant self-absorption of the scintillation light. BGO can be machined to various shapes and geometries. The crystal housing can be simple since no hermetic sealing is required.

The scintillation emission maximum of BGO is situated at 480nm. Figure 1 shows the emission spectrum. The light emission in photons/keV is about 15-20% of NaI(Tl); but, since the emission is partly in the area above 500nm where phototubes are less sensitive, the relative photoelectron yield of a bialkali PMT compared to NaI(Tl) amounts to 10-15%.

Figure 2 shows a pulse height spectrum obtained by irradiating a BGO crystal with 662 keV γ -rays.

Due to the high Z value of the material, the photofraction for γ -ray absorption

is high; and BGO scintillation crystals are used in applications where a high photofraction is required (for example, PET scanners) or because of its high detection efficiency (for example, Compton suppression spectrometers). It is a combination of properties that make BGO the material of choice for neutron activation analysis. Figure 4 shows the photopeak efficiency (also called the photofraction) – the ratio of the number of counts in the total absorption photopeak to the total number of counts as a function of the γ -ray energy for 38mm diameter, 38mm high (1.5" x 1.5") NaI(Tl) and BGO crystals.

The decay time of BGO is about 300ns at room temperature, which is comparable to that of NaI(Tl). As there is no slow component in BGO and the rise time is quite fast (intrinsic scintillator), it is possible to get good timing <2ns with 3" thick crystals .

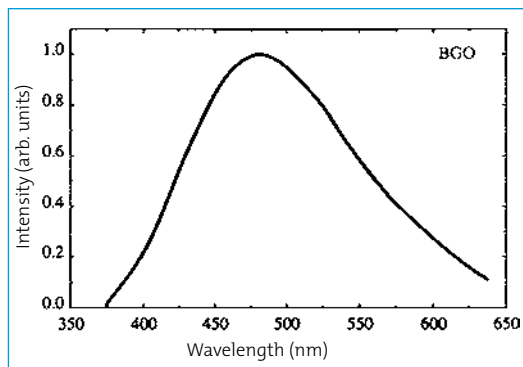


Figure 1. Scintillation emission spectrum of BGO

Properties

| | |
|---|----------------------|
| Density [g/cm ³] | 7.13 |
| Melting point [K] | 1323 |
| Thermal expansion coefficient [C ⁻¹] | 7 x 10 ⁻⁶ |
| Cleavage plane | none |
| Hardness (Mho) | 5 |
| Hygroscopic | no |
| Wavelength of emission max [nm] | 480 |
| Lower wavelength cutoff [nm] | 320 |
| Refractive index @ emission max. | 2.15 |
| Primary decay time [ns] | 300 |
| Light yield [photons/keV γ] | 8 - 10 |
| Photoelectron yield [% of NaI(Tl)] (for γ -rays) | 15 - 20 |
| Temperature response | -1.2%/C |
| Neutron capture cross-section | 1.47b |
| Afterglow @ 20ms | 150ppm |



BGO Bismuth Germanate Scintillation Material

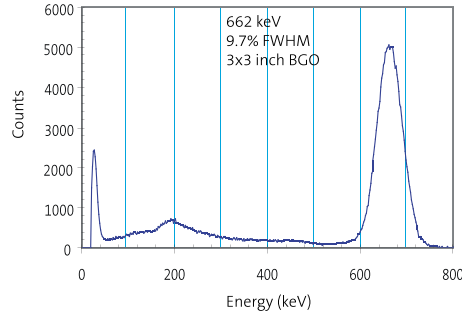


Figure 2. The pulse height spectrum of a 3" diameter, 3" high BGO crystal when irradiated with 662 keV γ -rays

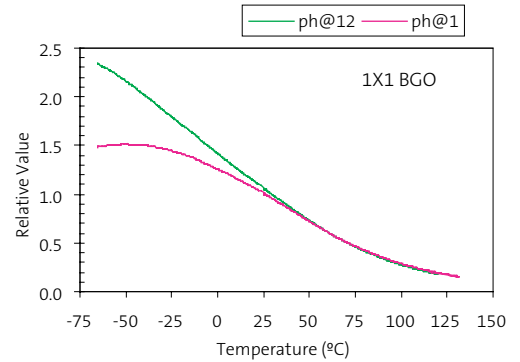


Figure 3. Temperature dependence of the scintillation light output of BGO (1 μ s and 12 μ s are the main amplifier shaping times)

The scintillation intensity of BGO is a strong function of the temperature. Figure 3 shows the relation. At room temperature, the rate of change with temperature is approximately -1.2%/C.

The radioactivity in BGO can make BGO unacceptable for some applications. We have developed a production process that significantly reduces the natural background, making our BGO well-suited for most applications.

BGO scintillation crystals are susceptible to radiation damage starting at radiation doses between 1 and 10 Gray ($10^2 - 10^3$ rad). The effect is largely reversible with time or annealing. Since the radiation damage to BGO crystals depends on the presence of sub ppm impurities, large differences between individual crystals can occur.

It is possible to read out BGO crystals with silicon photodiodes but, due to the moderate light output, this is only useful for the detection of high energy particles, or photons of more than a few MeV.

USA

Saint-Gobain Crystals
17900 Great Lakes Parkway
Hiram, OH 44234
Tel: (440) 834-5600
Fax: (440) 834-7680

Europe

Saint-Gobain Crystals
104 Route de Larchant
BP 521
77794 Nemours Cedex, France
Tel: 33 (1) 64 45 10 10
Fax: 33 (1) 64 45 10 01

P.O. Box 3093
3760 DB Soest
The Netherlands
Tel: 31 35 60 29 700
Fax: 31 35 60 29 214

Japan

Saint-Gobain KK, Crystals Division
3-7, Kojimachi, Chiyoda-ku,
Tokyo 102-0083 Japan
Tel: 81 (0) 3 3263 0559
Fax: 81 (0) 3 5212 2196

China

Saint-Gobain (China) Investment Co., Ltd.
15-01 CITIC Building
19 Jianguomenwai Ave.
Beijing 100004 China
Tel: 86 (0) 10 6513 0311
Fax: 86 (0) 10 6512 9843

www.detectors.saint-gobain.com

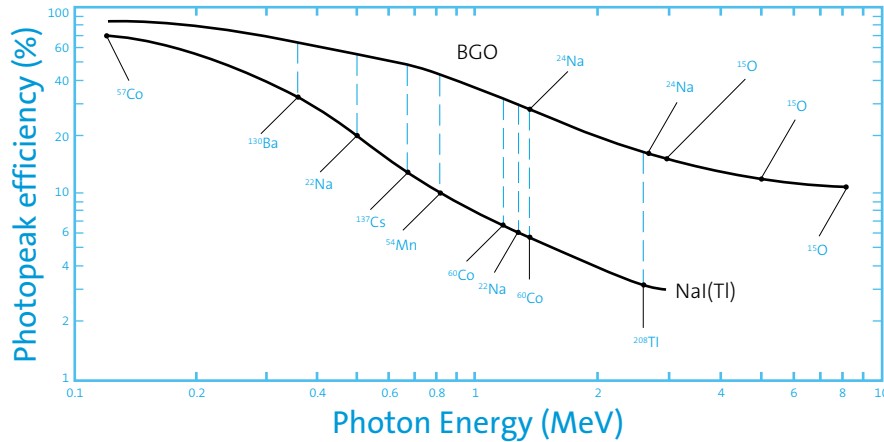


Figure 4. Photopeak efficiencies for BGO and NaI(Tl) scintillation detectors, 38mm diameter, 38 mm high.

Manufacturer reserves the right to alter specifications.

©2004-12 Saint-Gobain Ceramics & Plastics, Inc. All rights reserved.

(06-12)